

# IMPORTANT INFORMATION TO READ and RETURN

## Installation Requirements for a Whitley M155 HEPA Anaerobic Workstation

Thank you for choosing one of our products for your laboratory. To enable our engineers to perform an efficient, trouble-free installation please study, complete and email this form to us at [service@dwscientific.co.uk](mailto:service@dwscientific.co.uk). Should you have any questions, please do not hesitate to contact us, as we are here to help. When we have received the completed form, our Service Department will contact you to arrange a mutually convenient installation date.

**The following information represents the ideal requirement.  
Please contact us IMMEDIATELY if your intended location does not match this specification.**



### Access Requirements

Please note that this product is **wider than a standard door** so please ensure that you consider access carefully.

### Space Requirements

The weight of the equipment is 350kg. This workstation comes complete with a bespoke trolley. Please ensure that sufficient floor space is allocated. If you wish to locate the workstation on a bench, please contact us to discuss the implications.

#### External Dimensions\*

Width	Depth	Height	Height Including Trolley
mm	mm	mm	mm
2875	1056	1063	1911

\* Please Note:

If the **Refrigeration System** is fitted, please allow a clearance of 770mm behind the workstation. If the **Heat Removal System** is fitted, please allow a clearance of 600mm behind the workstation. Some of this clearance is necessary for air circulation.

### Gas Requirements

The incoming gas supplies must be terminated near the right hand side of the main chamber and fitted with leak-proof taps and pressure gauges.


The gas lines to which the equipment is attached are the responsibility of the user and should be constructed, tested and maintained to the standards specified within the British Compressed Gasses Association (BCGA) Code of Practice CP4 (or international equivalent). Gas lines previously used for flammable gases must be purged prior to re-use.

Regulators should be fitted in accordance with the information contained in the table below and the various pressures strictly adhered to.

Gas Type	Connection Details	Cylinder Regulator Required	Regulator Outlet Range	Flow Rate
Hydrogen (10%) / Nitrogen (90%)	¼" BSP male fitting or connection for 6mm Polyurethane tubing.  Hydrogen Regulator – Two Stage – order Code <b>A01745</b>	Two stage	4 - 6 bar (60-90 psi)	20 litres per minute (dynamic)
Nitrogen	¼" BSP male fitting or connection for 8mm Polyurethane tubing  Nitrogen Regulator – Two Stage – order Code <b>A01748</b>	Two Stage	4 - 6 bar (60-90 psi)	Minimum 150 litres per minute (dynamic)
Carbon Dioxide	¼" BSP male fitting or connection for 8mm Polyurethane tubing  CO2 Regulator – Two Stage – order Code <b>A01747</b>	Two Stage	4 - 6 bar (60-90 psi)	10 litres per minute (dynamic)

Air	¼" BSP male fitting or connection for 8mm Polyurethane tubing  Air Regulator – Two Stage – order Code <b>A01554</b>	Two Stage	4 - 6 bar (60-90 psi)	Minimum 20 litres per minute (dynamic)
-----	---	-----------	-----------------------	--

**Connection Type**



Push in connection

**OR**

Push on Connection

**Mains Requirements**

<b>Electricity Supply</b>	<b>Wall Socket</b>
240 volts	1 x Three Pin, 13 Amp. Minimum 6A rating

**Other Considerations**

Although the workstation should be located in a well ventilated area, avoid proximity to air conditioning systems and draughts caused by windows and doors.

Remember, if you do not have the required regulators you can order them from Don Whitley Scientific: :

- Hydrogen Regulator – Two Stage – order Code **A01745**
- Nitrogen Regulator – Two Stage – order Code **A01748**
- CO2 Regulator – Two Stage – order Code **A01747**
- Air Regulator – Two Stage – order Code **A01554**

**Notes**

If an existing unit is being taken in part exchange or is being removed from the laboratory, it must be de-contaminated before DWS staff handle the unit. A certificate or signed letter confirming the unit has been decontaminated must be given to our engineer.

In the UK, delivery and installation are free of charge (unless otherwise agreed). If our engineers are unable to install the unit and a return journey is necessary, **a charge may be made**. Export customers should please refer to their local distributor.

**It is essential that this form is completed and returned, to avoid delay to your installation.**

**THANK YOU FOR THINKING WHITLEY**

<b>Signature</b>	<b>Title</b>
<input style="width: 90%;" type="text"/>	<input style="width: 90%;" type="text"/>
<b>Print Name</b>	<b>Establishment</b>
<input style="width: 90%;" type="text"/>	<input style="width: 90%;" type="text"/>